

MCA8000D Multi Channel Analyzer

Features

- · Compatible with traditional analog pulse shaping
- MCA and MCS modes
- High speed ADC (100 MHz, 16 bit) with digital pulse height measurement
- 8k data channels
- · Minimum pulse peaking time 500 ns
- Conversion time 10 ns
- Sliding-scale linearization
- Differential nonlinearity <±0.6%
- Integral nonlinearity <±0.02%
- Two peak detection modes: first peak after the threshold (nuclear spectroscopy) or absolute peak after the threshold (particle counter calibration in clean rooms)
- Two TTL compatible gates for coincidence and anticoincidence
- Supports USB, RS-232, and Ethernet communication interfaces
- Dimensions: 5 x 2.8 x 0.8 in / 125 x 71 x 20 mm
- Weight: <165 gUSB powered
- Free Display Acquisition Software supports ROI, energy

- calibration, peak information, MCA configuration, and file management
- Free Software Developer's Kit (SDK)
 Complete protocol and ex- ample code to be used in custom software applications
- Optional PA Package for airborne and suspended particle detection applications



Amptek's MCA8000D is a state-of-the-art, compact, high performance, digital MultiChannel Analyzer (MCA). It replaces the MCA8000A, provid- ing the same functionality with reduced dead time, modern communications interfaces (USB, RS-232, and Ethernet), and much higher data transfer rates. Because it does not implement digital pulse shaping, it is compatible with analog pulse shaping systems.

Superior performance is obtained from a high speed ADC (a 100 MHz 16 bit ADC) and sophisticated digital logic. The MCA8000D digitizes the input signal (the output of an analog shaping amplifier) to measure the pulse height and obtain the pulse height spectrum. The high speed ADC, with a 10 ns conversion time, reduces dead time thus increasing throughput. The digital circuitry yields high accuracy and stability.


The MCA8000D includes three interfaces, USB (USB 2.0 full speed), RS-232 (up to 115 kB), and Ethernet (10base-T). It uses the same DPPMCA software package as Amptek's digital signal processors, including the DP5, PX5, and X-123. The MCA8000D can be powered over USB or with an external AC/DC power supply.

Compromising nothing in performance, the MCA8000D is a low power, light weight instrument which is exceptionally versatile and easy to use.

Fits in a shirt pocket!!

The **MCA8000D** is a full-featured digital multi- channel analyzer intended to be used with a wide variety of detector systems.


Examples of unipolar (left) and bipolar (right) input pulse shapes suitable for the MCA8000D

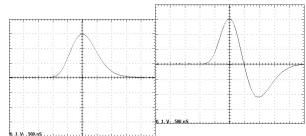


METEK

Wor.	
wer	
perating Power 2 W.	
wer Sources Power: +5V at 0.4A	
Input range: +4V to +5.5V (0.5 to 0.4A	
typical) Power source: USB power or	
AC/DC adapter (supplied)	
nnections	
The analog input accepts positive unipolar	•
or bipolar semigaussian type pulses of shaping time constants ≥200 ns or peakin	~
time	y
≥ 500 ns. See Figures 1 and 2.	
The dynamic range is 0 to +1 V or 0 to +10)
V, software selectable. Minimum pulse	-
height is 5 mV.	
Input impedance is 1 kΩ or 100 kΩ	
· · ·	
(options at purchase). The input has	
overload protection up to ±20 V.	
The DC level of the input signal must be	
tes The MCA-8000D has two logic gates. GATE	1
is generally used to synchronize acquisition	
	is
generally used for pileup rejection.	
USB: Standard USB Mini-B jack. Ethernet:	
rerface (I/O) Standard Ethernet jack.	
RS232: Standard 2.5 mm stereo audio	
jack. Power Center positive 5.5 mm x 2.1	
mm plug.	
nanical	
eight <165 g	
mensions 5 x 2.8 x 0.8 in (125 x 71 x 20 mm).	
formance Characteristics	
High speed 100 MHz, 16 bit ADC with	
Ise-Height digital pulse height measurement.	
Input ranges of 0 to 1V or 0 to 10V	
(software selectable) Number of channels	
commandable to 256, 512, 1k, 2k, 4k, or 8l	<
channels	

Pulse Processing	Pulse peaking time ≥ 500 ns to meet
Times	specifications.
Peak Detection	First Peak After Threshold (V1)
Modes (See	, ,
Figure 3)	
	Absolute Peak Above Threshold (V2)
Acquisition Modes	MCA and MCS (multichannel scaling) down
	to 10 ms/channel.
Differential	$<\pm0.6\%$ from 15 mV to full scale.
Nonlinearity	
Integral	<±0.02% over full scale.
Nonlinearity	
Gain Stability	Gain ±10 ppm/°C.; Zero drift ±10 ppm/°C. (typical)
Low Level	Software selectable threshold, in
Discriminator	increments of one channel
Counts per	Maximum = 16.7×10^6 counts (3 bytes).
Channel	iviaximum = 10.7 x 10 counts (3 bytes).
Real and Live	Data acquisition real or live time preset in multiples of one second. Maximum preset time 16777215 (2 ²⁴ - 1) seconds. Shorter or
Timers	non-integer accumulation times can be achieved by gating the input signal with Gate 1 or Gate 2.
	Real and Live Time: 4,294,967.29 s (0.01 s precision); 999,999.999 s (0.001 s precision). Acquisition Time: 99,999,999.9 s (0.1 s precision) Preset Counts: 4,294,967,295 (2 ³² -1) in channel or ROI
	RS232: 115 Kbaud.
Interface	USB: Standard 2.0 full speed (12
interrace	Mbps). Ethernet: Standard 10
	base-T.
Operating	-20°C to +60°C.
Temperature	
Exterior Controls and	Indicators
Status Light	Steady red light: power is ON and no data acquisition in progress. Flashing red light: data acquisition mode.

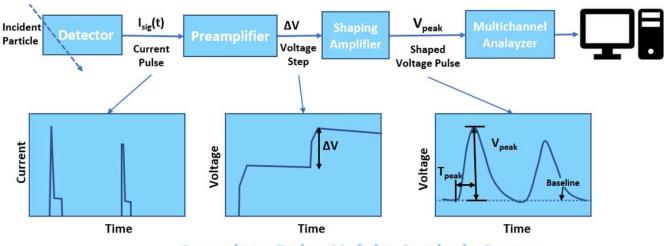

Complete Pulse Height Analysis System

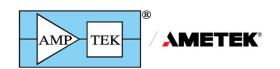

OPERATING NOTES

Amptek's MCA8000D is one component in a complete pulse height analysis system. Figure 3 illustrates a typical system: the radiation detector is connected to a charge sensitive preamplifier, which is connected to a shaping amplifier. The output of the shaping amplifier is a pulse with peak amplitude proportional to the energy deposited in the detector. The MCA measures the amplitude of the

peak and produces a histogram representing the pulse height spectrum. This histogram, an integer array of the number of events in each pulse height channel, is the primary output of the MCA and is transferred to a computer over USB, Ethernet, or RS232 interfaces.

Amptek's MCA8000D does not include pulse shaping circuitry but only the MCA circuitry. The MCA8000D must be preceded by pulse shaping circuits. Figures 1 and 2 show typical pulse shapes in- put to the MCA8000D. Amptek, Inc has other products (including the DP5, DP5G, PX5, and PX5-HPGe) which include both pulse shaping and MCA circuitry. These use digital pulse processing for the very best performance. Please contact Amptek, Inc for information on these inte- grated systems.



Note that some manufacturers use the term "MCA" for modules which include both the pulse shaping and the traditional MCA functions.

Amptek's MCA8000D is most often used to measure the energy spectrum from a radiation detector, where the pulse height is proportional to energy but can be used in other applications. For example, instruments used to monitor cleanrooms produce an output pulse proportional to the size of the particle, so the MCA8000D is used to measure the particle size spectrum. Some radiation detection systems used in timing produce a pulse height proportional to the time between events (using a time to amplitude converter), so the MCA8000D yields a timing spectrum.

SOFTWARE

The MCA8000D can be controlled by the Amptek DPPMCA display and acquisition software. This software sets the configuration (number of channels, input range, etc), can start and stop acquisitions, can display the spectra, and supports basic analysis functions such as defining regions of interest (ROI), performing calibrations, peak searching, and so on. Data are saved in an ASCII file format. The DPPMCA software is also used with Amptek's digital pulse processors, including the DP5, Gamma-Rad5, PX5, etc.

The MCA8000D comes with a complete Software Development Kit (SDK). The user can use this platform to easily develop software to control the MCA for custom applications or to interface it to a larger system. Examples are provided in VB, VC++, etc. The standard DPPMCA software runs only on Windows operating systems but users can develop interface software running on Linux or other operating systems using documentation and sample code which are provided.

ADVANCED ANALYSES

The pulse height spectrum is the primary output from the MCA8000D. It can also be used in a "multichannel scaling" mode to record count rate vs time (down to 10 ms/channel).

OPTION PA FOR PARTICLE COUNTING

The Option PA package has been developed to facilitate the use of the MCA8000D for particle counting in airborne (Size Calibration) and liquid suspended¹ (Number Calibration) particle applications. The unit is calibrated and certified traceable to the National Institute of Standards and Technology (NIST).

The Option PA package is capable of detecting pulses from 5 mV to 10 V. The MCA8000D is typically connected to the output of a particle sensor. It detects and displays a spectrum of pulse heights allowing the user to determine if a given particle size is producing the correct voltage.

The software included with the MCA provides information on the peak center (centroid and mean calculation) making it easy to determine if the peak is in the correct position. The supplied calibration curves convert the MCA channel scale to a mV scale. The calibration curves can be set to load automatically when the software opens.

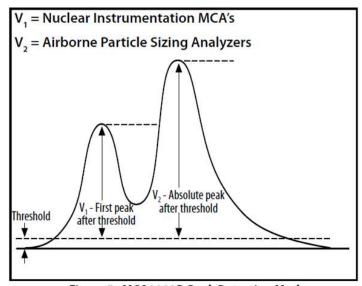


Figure 5. MCA8000D Peak Detection Modes

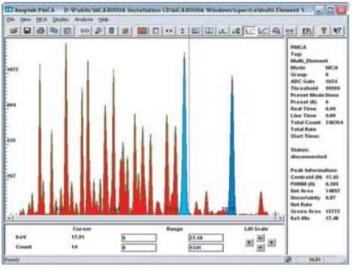


Figure 6. Example of DPPMCA Software for MCA8000D

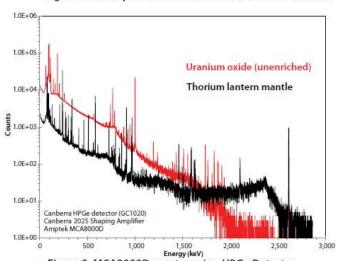


Figure 8. MCA8000D spectra using HPGe Detector

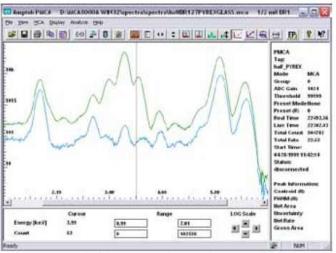


Figure 7. Example of DPPMCA Software showing multiple spectra capability

Frequently Asked Questions

How does the MCA8000D differ from the previous MCA8000A?

- 1) It now supports USB, Ethernet, and RS232 interfaces with much faster data transfers. It makes operation with modern computers very easy.
- 2) It uses the DPPMCA acquisition software, a new version of the older ADMCA software.
- 3) It can be powered by USB or an external plug. It does not contain a battery.
- 4) It has 8k data channels and is specified for pulse risetimes >0.5 us.
- 5) It is an all digital solution, with a fast ADC and digital peak detection.
- 6) It is smaller and lighter.


MCA8000D Ordering Information

MCA8000D Basic Package - for **Nuclear Instrumentation**

- MCA8000D
- One AC adapter for MCA8000D
- One BNC signal cable (1 meter length)
- One Mini-USB cable
- DPPMCA Display Software installation disk

MCA8000D PA Package - for Airborne Particle Sizing

 MCA8000D Basic Package, modified for 100 $k\Omega$ input impedance

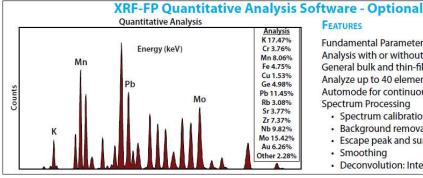

FEATURES

Fundamental Parameter (FP) Calculations Analysis with or without standards General bulk and thin-film analysis Analyze up to 40 elements Automode for continuous or repeated analyses Spectrum Processing

- Spectrum calibration
- Background removal and blank subtraction
- · Escape peak and sum peak removal
- Smoothing
- · Deconvolution: Intensity extraction



Certificate of Calibration



14 DeAngelo Drive, Bedford, MA 01730-2204 USA +1 781 275-2242 *Fax:* +1 781 275-3470 sales@amptek.com www.amptek.com